# Differential Plane-Mirror Interferometer



# **SP-DI Series**



### **Design and Operation**

SP-DI Series diifferential interferometers are used for highly precise differential length and angle measurement.

The strictly symmetrical optical structure of the interferometer achieves extremely high long-term stability of the length measurement.

Two parallel beams detect the relative motion between a reference point and the measuring point with the highest resolution and precision. The beam distance calibrated in the factory enables angles to be measured with high precision.


The measured values are recorded and displayed on a PC with optional data acquisition and display software.



#### **Major Performance Features**

- Differential length and angle measurement with the highest accuracy
- · Sensor head is made of stainless steel as standard
- · High long-term stability
- Differential measurement minimizes environmental effects
- · Easy adjustment and handling
- HeNe laser with high frequency stability as the metrological standard
- · Fiber-optic coupling of the sensor head
- Correction of environmental influences on the wavelength of the laser light
- Open interfaces for OEM software under Windows and Linux
- Other beam distances are possible

#### **Operating Principle**

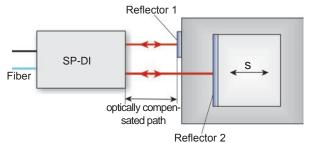

## **Applications**

- Highly precise differential length measurements, for example on positioning systems, for long-term material investigations and in dilatometry
- · Angles and tilts are measured with the greatest accuracy by calibrating the beam distance
- · Optional vacuum-compatible designs

Technical Data		Model SP 120 DI*	Model SP 2000 DI	
Measurement range	mm	70	2000	
Resolution	pm	20	20	
Laser waverlength	nm	632.8	632.8	
Frequency stability of the HeNe laser (after warm-up time)		≤ 3 · 10 <sup>-7</sup>	≤ 2 · 10 <sup>-8</sup>	
Warm-up time of the HeNe laser	min	1	1020	
Beam distance (standard)	mm	21		
Angular measurement range	arcmin	±1.5		
Angular resolution at 0.1 nm length resolution	arcsec	0.001		
Operating temperature range	°C	1530		
Maximum displacement speed of the measuring reflector	mm/s	800		
Dimensions (L x W x H)				
Sensor head with base plate	mm	180 x 140 x 43		
Electronic supply and evaluation unit	mm	450 x 400 x 150		
Mass				
Sensor head with base plate	kg	3.3		
Electronic supply and evaluation unit	kg	ca 8		
Interfaces standard		RS232C, USB		
optional		Digital 32-bit pa	Digital 32-bit parallel interface	
		Digital incremental signals (TTL level)		
		Analog incremental signals (1V <sub>PP</sub> )		
Cable length between sensor head and electronics unit	m	3, optionally up to 10		
Line voltage / frequency	VAC/Hz	100240 /4760		
Laser safety class according to EN 60825-1:2007 and		2M		
ANSI Z136.1 (CDRH)		I	I	

\*For ultra-stable measurements we recommend the model SP 2000 DI due to higher frequency stability of the HeNe laser.

# SIOS Meßtechnik GmbH

Am Vogelherd 46

98693 Ilmenau, GERMANY

Warning:

